

UltraTag™ Multiplex Labeling Kit Protocol — Chromogenic

The UltraTag Multiplex Labeling Kits efficiently conjugate 50 µg of antibody in 50 µL with one of three available Tags in two hours, with highly efficient tagging, and >80% recovery. Antibodies labeled with different UltraTag kits (UTCK-HRP1, UTCK-HRP2, UTCK-AP1) can be combined to create three-plex chromogenic multiplex IHC panels using included enzyme-labeled anti-Tag antibodies.

The UltraTag Labeling Kit modification protocol is summarized below:

1. Buffer exchange 50 µg carrier-free antibody into Mod Buffer.
2. Add buffer-exchanged antibody to Antibody Linker.
3. Add linker-modified antibody to UltraTag Reagent tube.

The only equipment required are a 200 µL pipette and a microcentrifuge.

This protocol can be downloaded from cellidx.com/documents/protocols/UltraTag-Multiplex-Labeling-Kit-Protocol-Chromogenic.pdf

Before Starting

- * Ensure your antibodies for labeling are at 0.9–1.1mg/mL in buffer free of extraneous proteins such as BSA, serum, or gelatin. Antibody formulations containing azide, trehalose, or other sugars, as well as any buffer containing tris, are acceptable.
- * Select antibodies that work well in your desired application. Antibodies with strong binding profiles are usually good candidates. We recommend selecting clones which stain strongly at less than 1 µg/mL when tested by conventional IHC. However, each antibody is different, and performance of individual clones cannot be guaranteed.
- * Enzyme-labeled rabbit anti-Tag antibody sufficient to stain 15–20 slides is included. Additional enzyme-labeled anti-Tag antibody is available. Components needed, but not supplied, include chromogens and HRP-arrest solution.
- * Optimization recommendations: For chromogenic multiplex staining, we recommend planning out the desired chromogen color for each antibody and tagging accordingly. For example, if you plan on using a Red AP chromogen for a specific antibody, we recommend tagging that antibody with UT015 since this kit comes with an anti-UT015 AP antibody.

Kit Contents

Section	Row	Description
A	1	BLUE -capped desalting column to buffer exchange antibody with collection tube
	2	Antibody-linker reagent for modification of 50 µg IgG @ 0.9–1.1 mg/mL
	3	GREEN -capped desalting column to buffer exchange linker-modified antibody with collection tube
	4	UltraTag Reagent
B	1	Link buffer
	2	RED -capped desalting column to buffer exchange Tag-antibody conjugate into PBS with collection tube
	3	Enzyme-labeled rabbit anti-Tag detector antibody, 15 µg @ 1.0 mg/mL
	4	Intentionally left empty



Antibody Preparation & Modification

1. **Desalting Column Preparation:** Break off bottom of **BLUE**-capped column, loosen the **BLUE** cap, and place column in collection tube, then place in microcentrifuge and spin for 2 minutes @ 1500 g. **Discard** flow-through.
2. Add your antibody solution (50 µg @ 1 mg/mL) directly to the **BLUE**-capped column, replace **BLUE** cap loosely on column, insert in a collection tube, and centrifuge for 2 minutes @ 1500 g.
3. Add flow-through from **BLUE**-capped column to Antibody Linker Reagent, pipette several times to mix, lightly vortex, and incubate for 1 hour @ room temperature.
4. **Desalting Column Preparation:** Break off bottom of **GREEN**-capped column, loosen the **GREEN** cap, and place column in a collection tube, then place in microcentrifuge and spin for 2 minutes @ 1500 g.
5. Add linker-modified antibody solution from Step 3 to **GREEN**-capped column, replace **GREEN** cap loosely on column, insert in a collection tube, and centrifuge for 2 minutes @ 1500 g.
6. Add flow-through from **GREEN**-capped column collection tube from Step 5 to UltraTag reagent tube, pipette several times to mix, and lightly vortex. Immediately continue to Step 7.

UltraTag Labeling & Purification

7. Add 7.5 µL of Link Buffer directly to the UltraTag reagent tube from Step 6, pipette several times to mix, lightly vortex, and incubate for 1 hour @ room temperature.
8. **UltraTag-conjugated Antibody Purification:** Break off bottom of **RED**-capped column and loosen the **RED** cap, place column in a collection tube, then place in microcentrifuge and spin for 2 minutes @ 1500 g. **Discard** flow through.
9. Transfer the UltraTag-antibody conjugation mixture from Step 7 to the **RED**-capped column and centrifuge for 2 minutes @ 1500 g.
10. Transfer the conjugate to a final storage tube of your choice. The conjugate yield of this process is ~70–80%. Therefore, the final concentration of antibody will be ~0.7–0.8 mg/mL.

Staining

- * Recommended concentration for staining with UltraTag-labeled primary antibodies is 5–10 µg/mL, however, optimization may be necessary, and performing a titration may suggest a different optimal concentration.
- * Recommended concentration for staining with enzyme-labeled anti-Tag detector antibody is 5 µg/mL.

The standard protocol for antigen retrieval and staining is comprised of the following steps:

1. **Antigen Retrieval:** Use your preferred antigen retrieval method for primary antibody/anti-species secondary antibody staining. See suggested protocols below

Antigen Retrieval Method	Buffer Recommended	Antigen Retrieval Time
Manual: Pressure Cooker	Citrate pH 6	15 minutes
Manual: PT Module	Dewax and HIER Buffer L (TA-100-DHBL)	5 minutes
BOND RX	ER2	20 minutes

2. **Staining:** Download detailed staining protocol from cellidx.com/documents/protocols/UltraTag-Multiplex-Staining-Protocol-Chromogenic.pdf

Detailed protocols for manual antigen retrieval and fluorescent multiplex IHC staining can be downloaded from cellidx.com/technical/protocols.



Imaging

The table below presents the Cell Palette chromogen catalog numbers and mounting media compatibilities of the chromogens which can be used with the enzyme-labeled anti-Tag antibodies included with the respective UltraTag kits. We recommend using CPAR-050 Cell Palette AP Red Chromogen, CPHB-050 Cell Palette HRP Blue Chromogen, and CPHY-050 Cell Palette HRP Yellow Chromogen. A combination kit (cat# CPK-015) of all three chromogens is available. For this chromogen combination, we have determined the optimal order to be Red, Blue, then Yellow. We recommend the use of a purple hematoxylin counterstain (cat# HP-050) followed by aqueous mounting medium or Cell IDx ChromoSealant (cat# CS-015) to seal stain followed by resin-based mounting medium.

Anti-Tag Antibody	Enzyme	Chromogen Recommendation	Catalog Number	Recommended Mounting Media
UT015	Alk Phos	Cell Palette AP Red	CPAR-050	Xylene-Based or Aqueous
UT016 or UT021	HRP	Cell Palette DAB	CPHD-050	Xylene-Based or Aqueous
UT016 or UT021	HRP	Cell Palette Blue	CPHB-050	Aqueous
UT016 or UT021	HRP	Cell Palette Yellow	CPHY-050	Xylene-Based or Aqueous

Disclaimer

The products offered here are for research use only. Any commercial application will require a license from Cell IDx. Cell IDx UltraPlex technology is patented and has multiple patents pending. Please contact Cell IDx for information regarding licensing information. Information in this manual is subject to change without notice and does not constitute a commitment on the part of Cell IDx. It is supplied on an “as is” basis without any warranty of any kind, either explicit or implied. This documentation is proprietary information and protected by the copyright laws of the United States and international treaties. The manufacturer of this documentation is Cell IDx.

Safety Information

- * **For Research Use Only.** Not for diagnostic or therapeutic use.
- * **WARNING – CHEMICAL HAZARD.** Some chemicals used can be potentially hazardous, and can cause injury or illness.
- * Read and understand the Material Safety Data Sheets (MSDS) available at www.cellidx.com before you store, handle, or work with any chemicals or hazardous materials.
- * Minimize contact with and inhalation of chemicals. Wear appropriate personal protective equipment when handling chemicals (*e.g.*, safety glasses, gloves, or clothing). For additional safety guidelines, consult the MSDS.
- * Check regularly for chemical leaks or spills. If a leak or spill occurs, follow the manufacturer’s clean-up procedures as recommended in the MSDS.
- * Comply with all local, state/provincial, and/or national laws and regulations related to chemical storage, handling, and disposal.