



## Product Specifications

### UltraPlex HI-2B Fluorescent Multiplex IHC Panel – Human CD68, CD163, Pan-CK, PD-L1

Description: UltraPlex fluorescent multiplex IHC kit to detect human CD68, CD163, Pan-CK and PD-L1  
 Cat. No.: HI02B-005  
 Application: IHC-P. Multiplex fluorescent IHC staining of human FFPE tissue sections.  
 Product Overview: All components required to stain 5 multiplex slides plus a single negative control slide with a simple two step protocol.

1° Antibody	Clone	1° Tag	2° Antibody	2° Fluor	Absorbance (nm)	Emission (nm)	Filter Set Examples	ExCoeff (M-1cm-1)
CD163	EPR19518	UT015	Anti-UT015	CL-490	491	515	FITC/GFP	73,000
PD-L1	73-10	UT014	Anti-UT014	CL-550	550	575	Rhodamine/TRITC/Cy3	150,000
Pan-CK	AE1/AE3	UT016	Anti-UT016	CL-650	655	676	Cy5	250,000
CD68	C68/684	UT019	Anti-UT019	CL-594	594	615	Texas Red	92,000

Optimized for use on FFPE tissue sections.  
 Suitable for spectral imaging.  
 Additional negative control (secondary reagents alone) are available, see NEG01B  
 Depending on your imaging scanner or microscope you may want to use Product HI02A with 490, 550, 650 and 750nm fluors instead.

Components: Vial 1: Antibody Diluent Solution  
 PBS, 1% BSA, 0.2% Tween 20, 15mM sodium azide  
 Vial 2: Protein Block Solution  
 PBS, 3% normal rabbit serum, 0.1% TritonX, 15mM sodium azide  
 Vials 3-6: Primary Antibody-Tag  
 Each 10ul at 0.5ug/ul in PBS, 15mM sodium azide, suggested dilution 1/100  
 Vials 7-10: Secondary Antibody-Fluor  
 Each 12ul at 0.5ug/ul in PBS, 15mM sodium azide, suggested dilution 1/100

Storage: 2-8°C

Shelf-Life: 6 months

Protocols: Download and use the following suggested protocols:

	Manual	Bond Rx
Antigen Retrieval	<a href="#">LINK</a>	<a href="#">LINK</a>
Staining	<a href="#">LINK</a>	<a href="#">LINK</a>

Notes: Suggested Mounting Medium: Fluoroshield plus DAPI, ImmunoBioScience  
 DO NOT USE: Vectashield Mounting Reagent, Vector Labs

Disclaimer: *For In vitro Research Use Only. Not for diagnostic or therapeutic use. Suggested applications of our products are not recommendations to use our products in violation of any patent or as a license under any patent of Cell IDx, Inc. Product may not be resold or modified for resale without prior written approval of Cell IDx, Inc.*