



Product Specifications

UltraTag Multiplex Labeling & Detection Kit - UTK-550

- Description:** Kit to label 50 µg of customer provided, purified primary antibody with CH014 hapten, and CL550 fluor labeled anti-CH014 detector antibody to stain 15-20 slides by IHC
- Cat. No.:** UTK-550
- Application:** Multiplex immunoassays including fluorescent immunohistochemistry (IHC), immunocytochemistry (ICC) and ELISpot. It is recommended to select an antibody that works well in the intended application.
- primary antibody must be purified at a concentration of 0.9-1.1 mg/ml in buffer free of extraneous proteins such as serum, BSA or gelatin. Antibody formulations containing azide, trehalose or other sugars as well as any buffer containing tris are acceptable
 - For fluorescence based multiplex IHC we suggest pre-qualifying your primary antibody by conventional IHC. Antibodies which stain well below 1 µg/ml are good candidates for UltraTag. Each antibody is different and performance of individual clones cannot be guaranteed.
- Components:** All components required to label 50 µg of purified antibody with a recommended starting concentration 0.9-1.1 mg/ml in carrier-free buffer to Cell IDx hapten CH014. The kit includes 3 required spin purification columns, antibody linker and lyophilized hapten. Also included is CL550 fluor labeled rabbit anti-CH014 detector antibody to stain 15-20 slides – 30 µl of concentrated antibody at 0.5 µg/ul, in PBS, 15mM sodium azide, suggested dilution 1/100

Anti-Hapten Antibody	Fluor	Absorbance (nm)	Emission (nm)	Extinction Coefficient (M ⁻¹ cm ⁻¹)	Recommended Exposure Time (ms)
CH014	CL550	491	515	73,000	100-300

- Storage:** 2-8°C
- Shelf-Life:** Not determined.
- Protocol:** [LINK](#)
- Notes:**

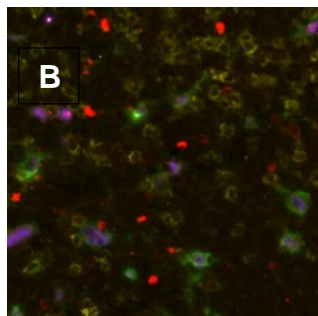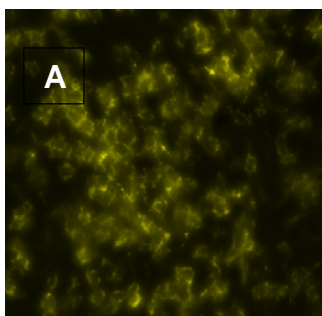

(A) CD4-CH014/anti-CH014-CL550, (B) UltraTag multiplex **CD163/CD4/CD68/LAG3**; primary and anti-hapten antibody staining @ 5 µg/mL; clones CD163 (EPR19518), CD4 (EP204), CD68 (68/684)/LAG3 (294)

Disclaimer: For in vitro Research Use Only. Not for diagnostic or therapeutic use. Suggested applications of our products are not recommendations to use our products in violation of any patent or as a license under any patent of Cell IDx, Inc. Product may not be resold or modified for resale without prior written approval of Cell IDx, Inc.